

Vista Quick Start Guide

A60-1017-551 ver. 1.0



A UNPACK YOUR BOX



1. Check your shipping carton and accessories case for the following items:
 - Vista instrument
 - Quick start guide
 - Power supply
 - Didymium diagnostic filter
 - HunterLab 2GB USB flash drive
 - Cleaning cloth
 - Stylus

Note: Please retain the packaging in case of instrument return to HunterLab.

2. Connect the power cord and all optional peripheral devices (mouse, keyboard, network cable, printer).



3. Turn the power on using the rocker switch on the right side of the instrument.



4. Once inside the software, the Color Data screen is automatically displayed.


B TAKING A MEASUREMENT

1. **STANDARDIZATION** – To begin, press the Workspace Icon in the upper right corner.

Workspace icon



2. Configure the Standardization Mode by selecting from the drop down menu. Check the "Include Haze" box if you plan to take both Color and Haze Measurements. Press "Standardize" to initiate this hands-free function.

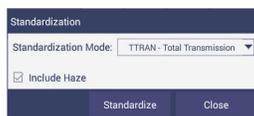

3. **READ** - Use the Read Measurement button to read samples. This button can be moved around by pressing and dragging to a new location.

Read Measurement icon



4. **VIEW FLIPPERS** --



These semi-transparent buttons are used to navigate between the measurement screens.

C CONFIGURE YOUR APPLICATION

To begin, press the WORKSPACE Icon in the upper right corner of the screen:

Workspace icon



1. **COLOR SCALES**-- Select and globally configure Color Scales/Indices /Differences.
2. **TOLERANCES**- Configure the Tolerances of the selected Scales, Indices and Differences.
3. **VIEWS**- Select the Color Measurement Screens.

D SAVE AND OUTPUT YOUR RESULTS IN "JOBS"

1. Press the JOBS Icon in the upper right corner.

Jobs icon



From the "JOBS" menu,

2. Select "Save Job" to store the your results in Vista's on-board memory.
3. Select "Data" to output results to a printer, email, or flash drive (CSV).

For more information, please refer to the manual.

E ACCESSORIES

A list of sample measurement accessories can be found at www.hunterlab.com/vista

F WARRANTY

HunterLab warrants that Vista is free from defects in material and workmanship under normal use and service. The time limit on this warranty is one year from date of shipment. For more information, see the Vista User's Manual.

For more detailed instructions on using the instrument, please refer to the Vista User's Manual.

Visit www.hunterlab.com for information on all our products.